

## 37000-006 Modular animal enclosure Assembly Instruction (Rev 1-Jan 2019)


To be used on Ugo Basile:

37450 Dynamic Plantar Aesthesiometer (DPA) and

37370 Plantar Test (Hargreaves Apparatus)

Your assembled enclosure will look like this:

A 37000-006, modular animal enclosure includes 3 modules, the parts for assembling the modules are in 3 separate boxes.

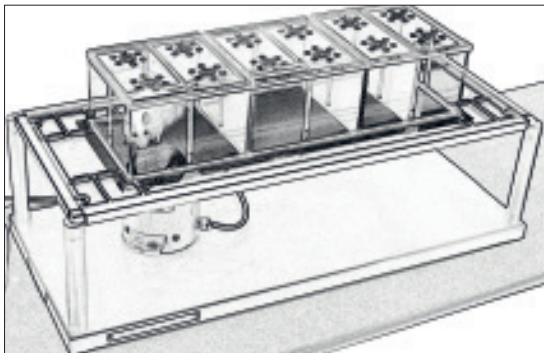

Whats in the box, for each cage module; quantity and codes:

**n.2 37000-006MD01**



**1**

**n.3 37000-006MD02**



**2**

**n.1 37000-006MD03**



**3**

**n.6 37000-006MD04**



**4**

**n.1 37000-006MD05**



**5**

**n.2 37000-006MD06**



**6**

**n.2 37000-006MD07**



**n.4 PLASTIC SPACERS**




## 37000-006 Modular animal enclosure Assembly Instruction (Rev 1-Jan 2019)

To be used on Ugo Basile:

37450 Dynamic Plantar Aesthesiometer (DPA) and

37370 Plantar Test (Hargreaves Apparatus)




How to assemble the enclosure:

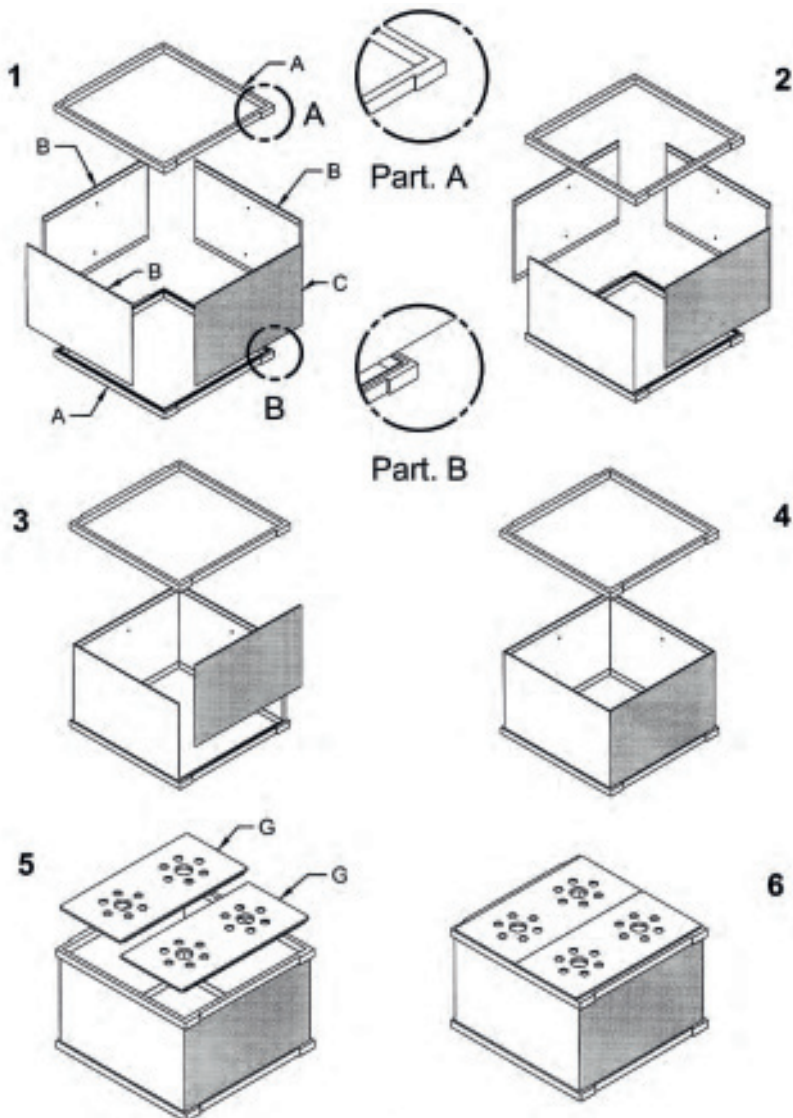


### 37000-006 Modular animal enclosure Assembly Instruction (Rev 1-Jan 2019)

To be used on Ugo Basile:  
37450 Dynamic Plantar Aesthesiometer (DPA) and  
37370 Plantar Test (Hargreaves Apparatus)


How to add the partitions:

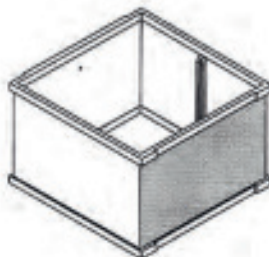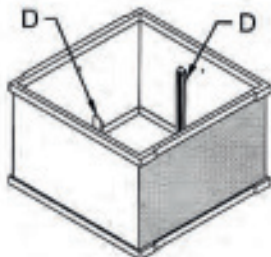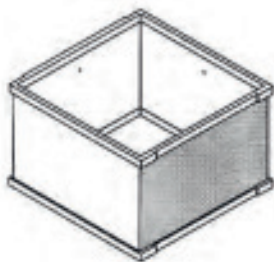


**4**

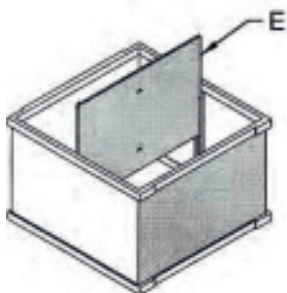


**5**

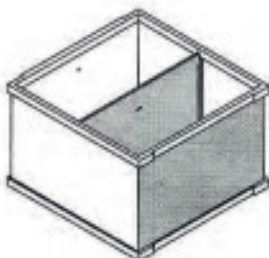


**6**

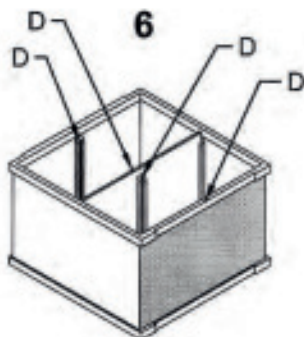


**7**

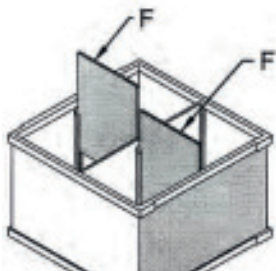

**8**

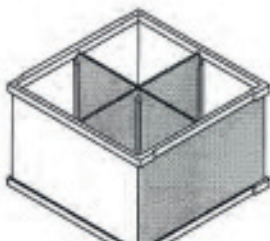

**9**

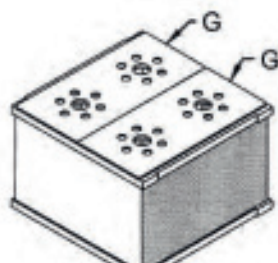



ugo basile®

TRANSFORMING IDEAS  
INTO INSTRUMENTS


## 37000-006 Modular animal enclosure Assembly Instruction (Rev 1-Jan 2019)

To be used on Ugo Basile:

37450 Dynamic Plantar Aesthesiometer (DPA) and

37370 Plantar Test (Hargreaves Apparatus)

How to assemble the single enclosure:

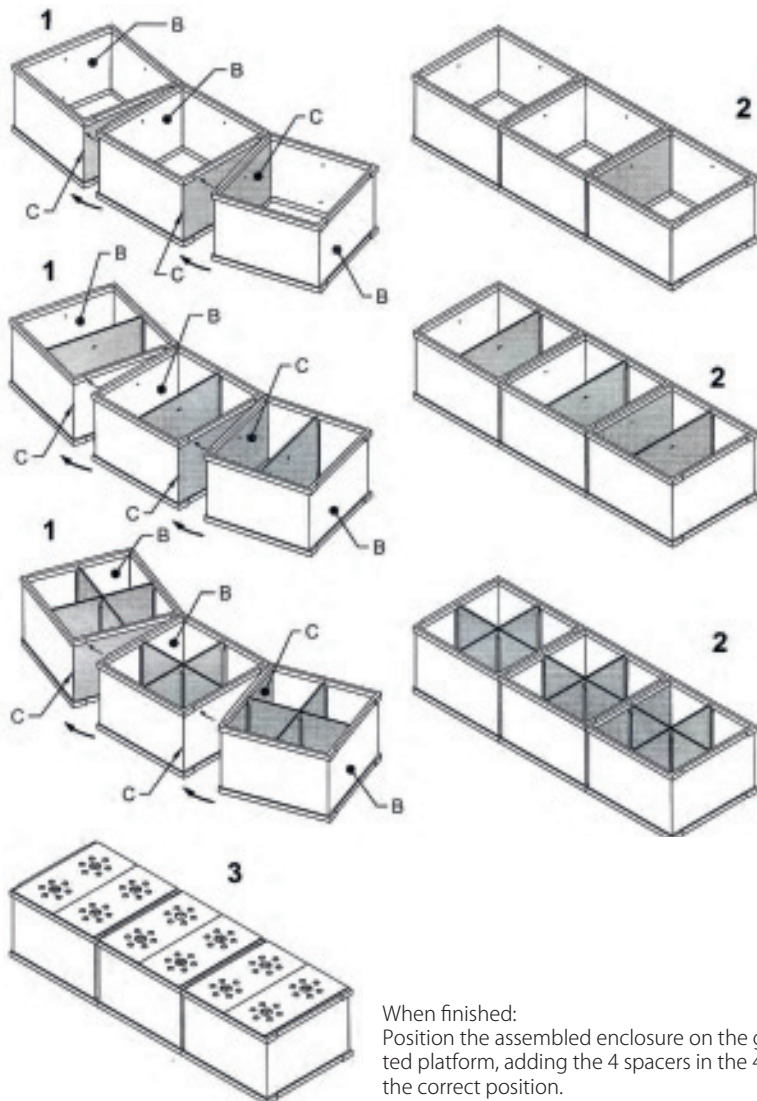

When finished:  
Position the assembled enclosure on the glass or on the perforated platform, adding the 4 spacers in the 4 corners, to keep it in the correct position.