


Mouse and Rat Restraints

Traditional Triangle Shape

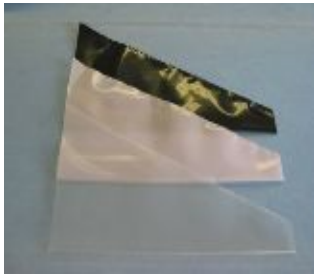


Includes instructions
& twist ties

Size	Animal	Clear	Black	Rose
X-Small <i>mice</i>	< 20 g	VI-1800C	xx	xx
Small	≤ 120 g	VI-1801C	VI-1801B	VI-1801R
Medium	≤ 225 g	VI-1802C	VI-1802B	VI-1802R
Large	≤ 450 g	VI-1803C	VI-1803B	VI-1803R

• 50 restraints/pack, custom sizes available •

Body-Contour™ Cone Shape

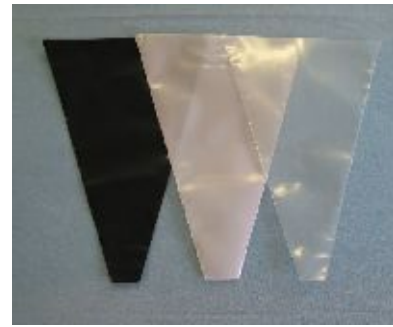

Size	Animal	Clear	Black	Rose
Small	≤ 120 g	VI-1801CV	VI-1801BV	VI-1801RV
Medium	≤ 225 g	VI-1802CV	VI-1802BV	VI-1802RV
Large	≤ 450 g	VI-1803CV	VI-1803BV	VI-1803RV

• 40 restraints/pack, custom sizes available •


Humane
Safe
Disposable

In Low Tail Injection Platform

Ref: VI-IL300



- ◆ Quick Restraint
- ◆ Low Stress
- ◆ Easy to Clean

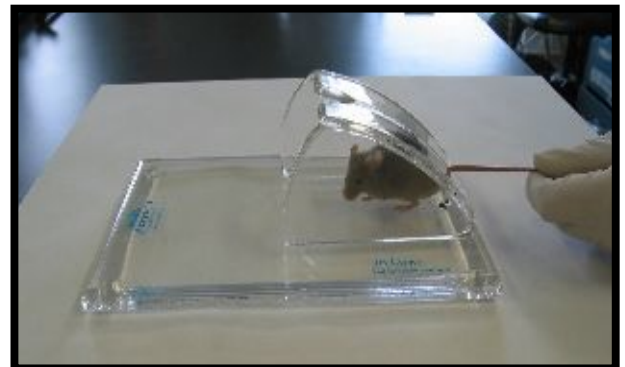

In Low Tail Cutting Platform

Cat #: VI-IL200



- X Quick Restraint
- X Tube Holder
- X Ruler for Accurate Snips
- X Easy to Clean



Rodent Restraint Bags




Step 1

Hold large end of bag open and allow rat to enter.



Step 2

As rat enters, keep narrow end of **bag & animal pointed up**.
Animals prefer to run upwards.

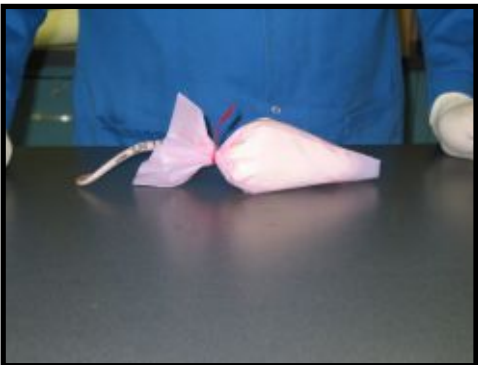

Step 3

Carefully twist plastic bag around base of tail and use twist tie to secure plastic to tail.

Technical Notes:

- Do not leave animal in bag for extended periods of time. Animal body heat will increase when restrained in a plastic bag.
- Placing a mark at the base of the tail (on top) before placing animal in bag will help with tail orientation when animal is fully immobilized.
- Black/Rose colored restraint bags reduce light and provide a more calming environment for animals.
- Body-Contour cone shape bags provide easier animal capture and increased animal comfort.